

Spectrafect® Diffuse Reflectance Coating

To maintain the unique optical and reflectance properties of Spectrafect Diffuse Reflectance Coating, the following handling procedures must be followed:

General Care and Cleaning

1. Wear powder free vinyl gloves whenever touching the coated surface.
2. Never attempt to clean the surface of a coated product with any form of liquid cleaner. Particles that contaminate the surface may be removed with a clean dry, filtered burst of compressed air. The air nozzle should be held no closer than 3 inches. Air compression should be no greater than 30 PSI.
3. Never place the coated surface face down onto another surface, unless it is specifically directed by a procedure.

For further information on the care and handling of Spectrafect coating, please contact the Labsphere Technical Sales Department.